


ELTRA ONH-2000 Simultaneous Oxygen / Nitrogen / Hydrogen Determinator

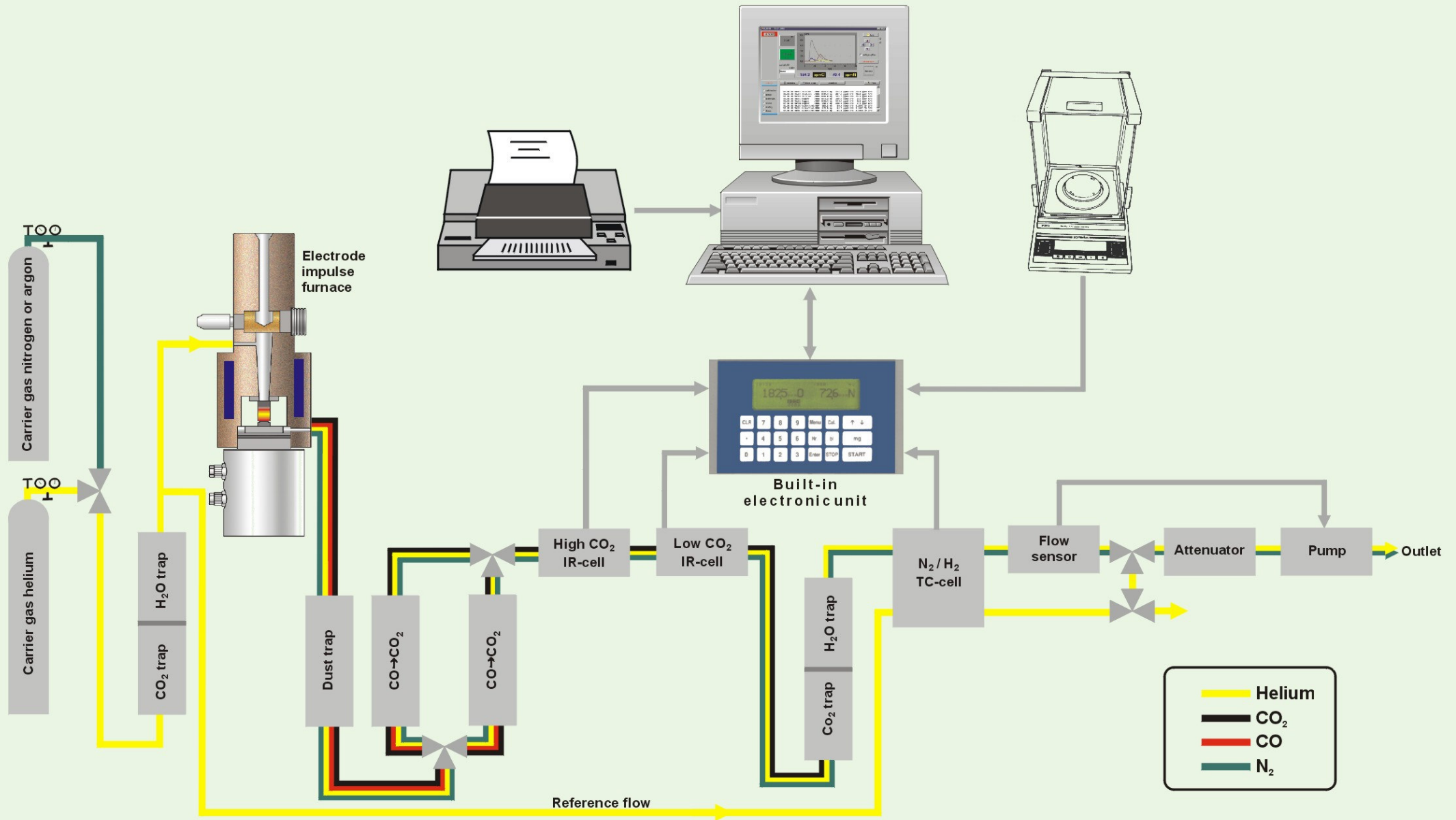

The ONH-2000 can alternatively operate as a simultaneous oxygen / nitrogen or as a simultaneous oxygen / hydrogen determinator. The changing from ON to OH mode and reverse, is easily done with a switch having an ON and an OH position. The carrier gas cylinders are also selected by solenoid valves controlled by the ON / OH switch. For OH operation, Schützes reagent is used as CO → CO₂ converter, while for ON operation the more economic CuO is used in a catalyst furnace.